



Klipspringer

Compliance with confidence

Rapid Allergen Test Kit INSTRUCTIONS


10 tests per pack




- ✓ Simple and easy-to-use (just 1-step).
- ✓ Specialized for swab and rinse water.
- ✓ No need for special equipment (Swab Kit is included)
- ✓ Limit of detection: 0.5ppm (μg protein/mL)
- ✓ Assay time: 10 minutes

Swab


1 Swab the specified area

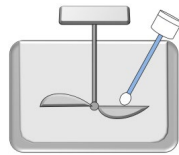

2 Shake vigorously

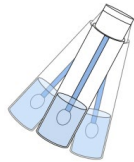


3 Add 200 μL (4 drops) to the test stick

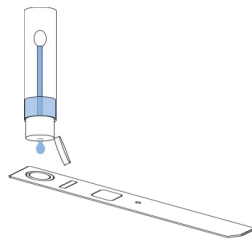


Rinse Water

1 Collect rinse water

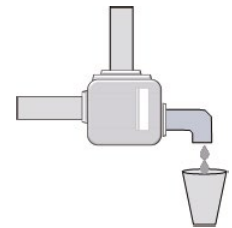

2 Add 200 μL (4 drops) to the test stick

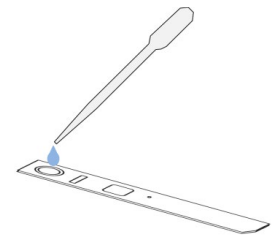

10 mins
to check the
result

Positive



Negative



Manufactured by

E-mail

Website

Morinaga Institute of Biological Science, Inc. (MIoBS), 2-1-16 Sachiura, Kanazawa-ku, Yokohama 236-0003, Japan

info_miobs_e@morinaga.co.jp

http://www.miobs-e.com

01473 461800

sales@klipspringer.com

www.klipspringer.com